

TT-4000 SMARTBOX® Tabletop Chamber

The TT-4000 Tabletop Chambers automate a three-stage euthanasia cycle at the push of a button: charging, dwell time, and exhaust. At the start of a cycle, the chamber lid automatically locks and will not unlock until CO₂ evacuation is complete. The clear acrylic construction provides full visibility and a top loading lid promotes easy cage access. The airtight, gasket-sealed lid prevents gas leakage during operation. The manifold design ensures even gas distribution. The system accommodates the full range of rodent cages from all manufacturers.

- **Compliant with AVMA Guidelines**
- Automation virtually eliminates operator error
- Lid automatically locks during operation
- After starting chamber, the technician is free to do other work
- Double throw switch allows for two different treatment times
- Programs can be customized based on facility requirements
- Operates at a fixed flow rate based on the external CO₂ source

Required Accessories (choose one):



EA-285
(use with cylinders)

or



1050-A1-BR
(use with piped-in gas)

Accessories



**EA-1040 Mobile
Battery Back-Up**




E-29000
(tanks not included)


SMARTBOX®




*TT-4000 includes:
SMARTBOX® Side Control Box, Tabletop Chamber,
Exhaust Blower, 8' of 4" Exhaust
Tubing, two 4" Hose Clamps
(Cart sold separately)*

Exhaust Blower included
Mounted when cart is purchased.

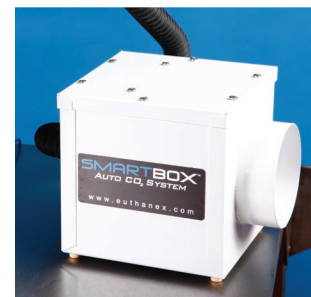


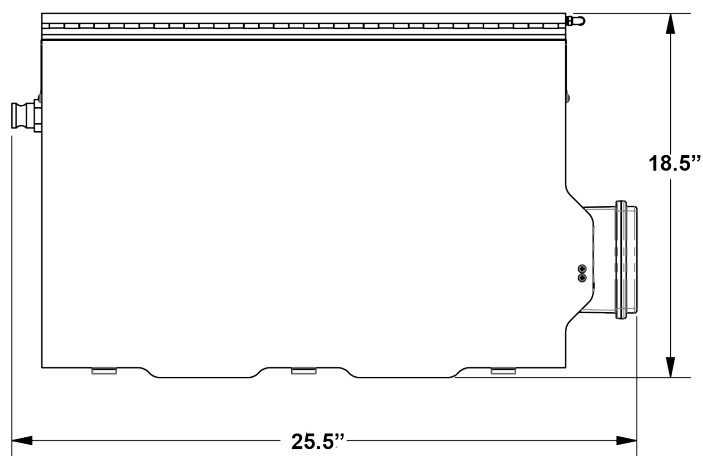


TOP VIEW

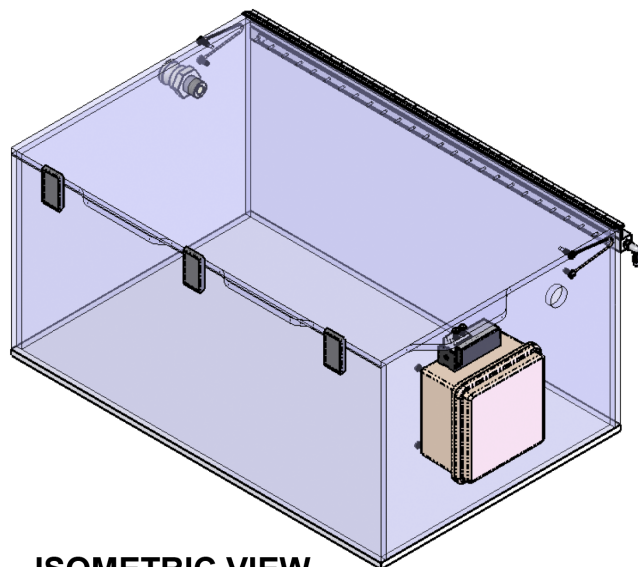

ISOMETRIC VIEW

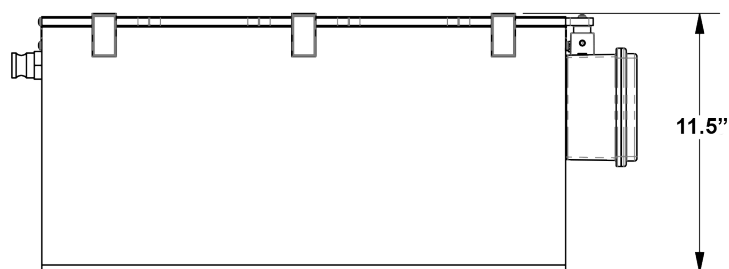

FRONT VIEW

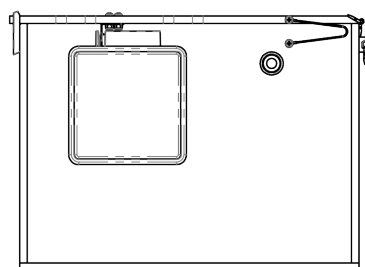

SIDE VIEW

INSIDE CHAMBER DIMENSIONS

20.75" L X 16" W X 10.5" H

REQUIREMENTS

ELECTRIC	110 VAC 15 AMP
FLOW RATE	19.0 LPM OR 40 CFH
4" DUCT	20CFM